User Manual 用户手册

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Thank you for using our products. As a company of caring employee safety and environmental protection, SYSBEL provides high quality products to ensure the safety of your working environment and personnel.

SYSBEL eyewash complies with the standard of ANSI Z358.1-2014. As the emergency safety protection equipment, eyewash station series can be used for flushing the parts exposed to the splashed chemicals and other substances, such as the body, face or eyes, to promptly reduce the damage caused by chemicals and achieve rapid primary self-aid.

It complies with the regulations in ANSI Z358.1-2014 standard and OSHA 29 CFR 1910.151(C) issued by ANSI has acquired CE Certification

Discription

WG6000A / WG6000B

- It adopts PE materials, and the production materials comply with the regulations of F.D.A on the grade on PE materials for food (F.D.A. Regulation on Polypropylene: the requirements of U.S.21 CFR F.D.A. Regulation Part 177.1520, portable containers clauses (C) for food contact use), and it is superior in its high intensity, good toughness, good Rigidity, heat and cold resistance
- It complies with human engineering principle in design, and adopts the law of gravity to produce water pressure and deliver water to two nozzles for eye washing
- It applies to work places without water supply and discharge facilities and can be installed and fixed onto walls as required to provide timely and professional protection
- It complies with regulations of ANSI/SEA on emergency eye wash station and showering equipment-ANSI Z358.1-2014
- It is suggested to discharge the liquid in eye wash station weekly or as required and clean every component of eye wash station
- Wall-mounted backing plate is equipped.
- Pull the rubber cover for immediate washing.
- Easily discharge water in tank with the drainage hole at the base.
- Two ABS spray heads with dust rubber covers.
- Three supporting feet make the eyewash steady and have enough space to operate when it is transported.

WG6000AD / WG6000BD

- Mobile waste cart adds portability and even holds the unit, also, collects waste water
- Perfectly equipped with WG6000A and WG6000B
- After placing on the cart, the spray heads are located between 83.8-134.6cm from the floor, which comply with ANSI Z358.1-2014

WG005

- Highly visible yellow, rugged polyethylene construction is well suited to long term use in harsh environments
- Built in handle for easy transport is great for use on remote site
- Petcock on the bottom for easy drainage
- Comes equipped with two rigid casters in front and two swivel casters with "Total Lock" brakes in back
- Includes clearly marked eyewash signs
- Cart has 37.5-gallon (142 liter) capacity and measures 86H x 61.5W x 52D (cm)

Trademark & Copyright

SYSBEL® is a registered trademark of SYSBEL, which is protected by law.

According to copyright law, it is forbidden to copy the whole or particle content of this manual without the written consent of SYSBEL. We have tried our best to ensure that the information in this manual is correct, and if the information is updated, we will not inform you. You can check our official website at www.sysbel.com.cn any time.

Portable Eyewash								
Product	Main Material	Water Capacity (L/Gal)	Flow Speed (L/min)	Flushing Time (min)	Packing size HxWxD(mm)	G.W. (kgs.)	Certification	Model
Type A	PE	60/16	2.6	18	436×750×435	9.5	ANSI Z358.1 AS 4775 CE	WG6000A
Type B		30/8	2.6	9	357×635×360	7		WG6000B

Portable Eye Wash Station							
Product	Water Capacity (L/Gal)	Flow Speed (L/min)	Flushing Time (min)	Matching Products	G.W. (kgs.)	Certification	Model
Type A	60/16	2.6	18	WG6000A+WG005	27.5	ANSI Z358.1 AS 4775 CE	WG6000AD
Type B	30/8	2.6	9	WG6000B+WG005	25		WG6000BD

Portable Eyewash with Cart						
Product	Main Material	Water Capacity (L/Gal)	Ext. size HxWxD(mm)	Packing size HxWxD(mm)	G.W. (kgs.)	Model
Mobile Waste Cart	PE	37.5/142	860×615×520	760×629×535	18	WG005

Instruction

Installation
STEP 1: According to ANSI Z358.1, make sure portable eyewash unit shall be installed in accessible locations requiring no more than 10 seconds to reach and shall be within less than 30 meters from the hazard. Employees shall be instructed in the proper use of emergency eyewash equipment.

Select a suitable mounting location for the WG6000A and the WG6000B capable of supporting 68kg and 38kg. Mount unit to wall (not designed for tabletop or shelf mounting) utilizes "J" bracket (supplied). Note: The 3/8" mounting hardware not included. The top of the mounting hole should be located between 143.8cm to 194.6cm from the floor. Mark the mounting locations onto the wall and place mounting bracket on surface. Drill necessary holes into mounting surface, and secure "J" brackets Mounting Bracket to surface using mounting hardware. The eyewash nozzles shall be between 83.8 cm (33 inches) and 134.6 cm (53 inches) above the floor when eyewash arm is in the active (lowered) position. (See the Structure Drawing).

IMPORTANT:
Filled eyewash unit (WG6000A) is heavy (approximately 150 pounds). Mounting bracket anchors, (Quantity-4, not supplied), or shelf must be strong enough to support the filled portable eyewash and the weight of a worker leaning on portable eyewash.

STEP 2: Equip with a waste water collection barrel, if desired.
STEP 3: Remove Fill Cap and fill the tank of portable eyewash with potable water (make sure that no debris or contaminants of any kind are present in it).
STEP 4: Take proper precautions to help prevent the growth of potentially harmful bacteria in eyewash tank. (Refer to Maintenance Instructions).

RECOMMENDED TOOLS: Four 3/8" diameter hardware suitable for supporting filled eyewash (approximately 150 pounds) and tools needed to install the hardware.

Operation

Instruct all employees in the use of this Emergency Eyewash. Personnel should be assigned to check eyewash liquid level daily.

STEP 1: Lead patient to nearest eyewash station. Seconds can make the difference in preventing serious eye injury.
STEP 2: Pull black eyewash arm firmly. Eyewash arm will drop to activate eyewash.
STEP 3: Lower head until eyes enter eyewash stream. Use thumbs and forefingers to hold eyelids open.
STEP 4: Seek qualified medical attention immediately

Maintenance

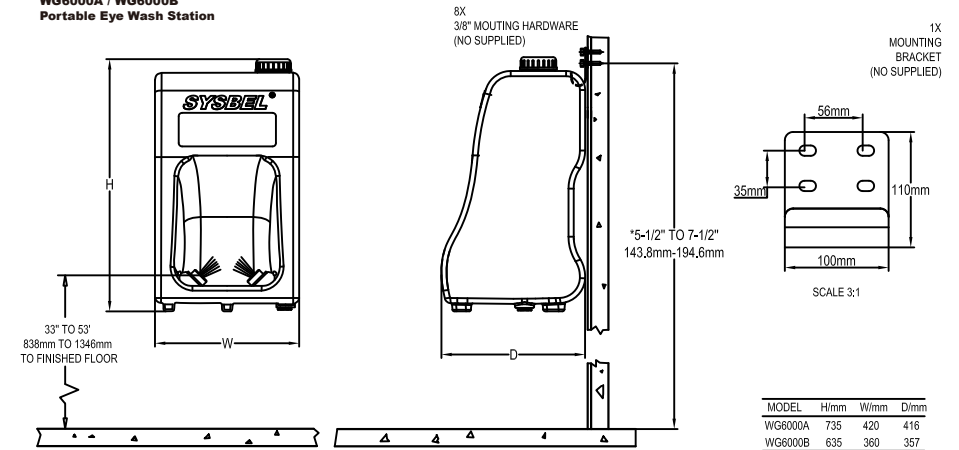
ANSI Z358.1 requires that all self-contained eyewash units shall be visually inspected weekly, unless the product is placed in environments with extreme conditions that may affect the form, fit, or function of the unit (temperature, dirt, etc.). If this is the case, then the inspection should be conducted more often as the end user deems necessary. Occasional cleaning and disinfection is recommended.

NOTE:


Customer must provide an inspection tag and maintain a record of inspection dates and type of maintenance performed
Take proper precautions to help prevent the growth of potentially harmful bacteria in eyewash tanks. We recommend the following procedure:
Drain, flush and refill portable units with clean potable water at least once a week. Thoroughly cleanse tank at least once every month.

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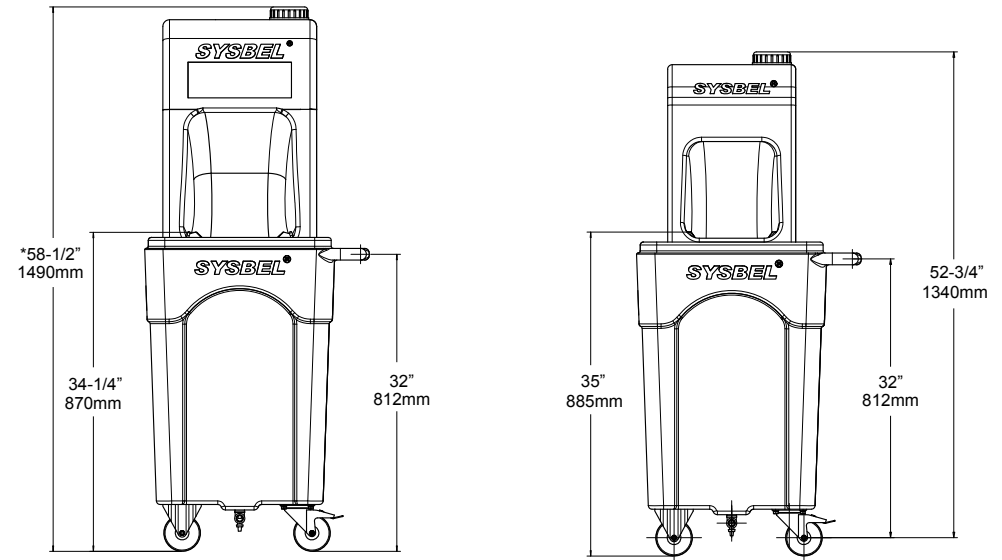NOTES:

- TO COMPLY WITH ANSI Z358.1-2014 AND OSHA 29 CFR 1910.151(C) FOR EMERGENCY EYEWASH:
 - UNIT SHALL BE INSTALLED IN ACCESSIBLE LOCATION REQUIRING NO MORE THAN 10s TO REACH AND SHALL BE WITHIN LESS THAN 30m FROM THE HAZARD.
 - EMPLOYEES SHALL BE INSTRUCTED IN THE PROPER USE OF EMERGENCY EYEWASH EQUIPMENT.
 - SELECT A SUITABLE WALL FOR THE WG6000Y CAPABLE OF SUPPORTING 150LB(68KG) TO LOCATE.
 - EYEWASH SHALL BE POSITIONED WITH NOZZLES NOT LESS THAN 838MM(33IN), AND ON MORE THAN 1346MM(53IN).
- UNIT MUST BE LEVEL FROM FRONT TO BACK AND SIDE TO SIDE.
- *DIMENSION MAY VARY 3/8"(9.525mm) ; TOLERANCE: ±8mm.